

Lovibond® Colour Measurement Tintometer® Group



Model F

Visual Colorimetry : Lovibond® Precision



- Glass Colour Standards Housed in Durable Racks
- Prismatic Optical System Provides Improved Colour Matching
- Balanced Field of View
- Replaceable Sample Chamber Liner

Part Number: 180000

Visual Colorimeter for Lovibond® RYBN Scale

The Model F is a visual colorimeter designed to optimise the use of Lovibond® glass filters in order to express colour in terms of Lovibond® Red, Yellow, Blue and Neutral.

Two Adjacent Fields of View

It is arranged with two adjacent fields of view, seen through the viewing tube, so that the product in the sample field and a white reflective surface in the comparison field are observed side by side, suitably illuminated.

Easy Colour Comparison

The Lovibond® glasses are introduced into the comparison field by a simple system of sliding racks, allowing you to compare the colour of light which is either transmitted through, or reflected from, the sample. The racks are varied until a visual colour match is found for the light from the sample and its colour can then be expressed in Lovibond® units.

Excellent Colour Matching

A series of neutral glasses in racks is also supplied; these can be introduced into the sample field to 'dull' the colour of products which are too bright to obtain a good colour match using Lovibond® Red, Yellow or Blue glasses. The racks allow a visual colour match to be found from the light from the sample and its colour can then be expressed in Lovibond® units.

Assured Accuracy of Lovibond® RYBN Colour

Lovibond® Tintometer® is the only company that can guarantee accurate, validated Lovibond® RYBN Colour readings.

Easy Maintenance for Optimum Performance

The Lovibond® Tintometer is a precision optical instrument which should be kept clean to preserve its accuracy, even in the harsh operating conditions often encountered. This is why the Model F has been designed to allow easy cleaning and maintenance with the minimum of cost and effort. The instrument includes a white plastic liner in the sample chamber which can be removed for cleaning if a spillage takes place. It can be replaced at intervals to maintain the whiteness of the interior, an essential feature for accurate colour matching.

Separate filters in the racks also simplify cleaning and allow replacement of single glasses (as opposed to complete racks) if required.

Application Versatility

The Model F is an easy-to-use instrument with versatile application in colour measurement of products which transmit light as well as opaque solids, powders and pastes (an optional 'solid' sample accessory pack may be required for colour grading of light-reflecting products).

Compliance to International Standards

It is available in slightly varying formats to meet the requirements of different applications and the national and international standardising bodies which specify the instrument in their official methods for colour measurement.

Industry

Chemical Industry | Food and Beverage Industry | Industries Others | Oil and Gas

Application

Beer, Malts and Caramels | Edible Oils and Fats | Food and Beverage | Petroleum Oils and Waxes

Model F

The Model F is a visual colorimeter designed to optimise the use of Lovibond® glass filters in order to express colour in terms of Lovibond® Red, Yellow, Blue and Neutral. It is available in slightly varying formats to meet the requirements of different applications and the national and international standardising bodies which specify the instrument in their official methods for colour measurement.

Technical Data

Applicable Standard	AOCS Cc 13e-92 AOCS Cc 13j-97 BS684 IS 548 (part 1)-2010 ISO 15305 ISO 27608 MS 252 : Part 16 MS 817 : Part 12
Color Type	Transparent Translucent Opaque
Path Length	Up to 153 mm (6")
Illuminants	Approximates to daylight
Portability	Benchtop
Compliance	CE
Dimensions	330 x 230 x 410 mm

Delivery Scope

- Model F Housing
- Complete set of 11 racks containing permanently coloured glass filters for the Lovibond® colour scale
 - (Red 0.1 - 0.9, 1.0 - 9.0, 10.0 - 70; Yellow 0.1 - 0.9, 1.0 - 9.0, 10.0 - 70; Blue 0.1 - 0.9, 1.0 - 9.0, 10.0 - 40; Neutral 0.1 - 0.9, 1.0 - 3.0)
- Sample chamber liner with a white PTFE powder reference
- Spare white reference
- A pair of spare bulbs
- 1 x 1" & 1 x 5¼" rectangular fused glass cells

Accessories

Title	Part Number
Lamp 12v10w	124100
Vacuum formed tray	180110
Halon Self Adhesive	183079

Tintometer GmbH
Lovibond® Water Testing
Schleefstraße 8-12
44287 Dortmund
Tel.: +49 (0)231/94510-0
sales@lovibond.com
www.lovibond.com
Germany

The Tintometer Limited
Lovibond House
Sun Rise Way
Amesbury, SP4 7GR
Tel.: +44 (0)1980 664800
Fax: +44 (0)1980 625412
sales@lovibond.uk
www.lovibond.com
UK

Tintometer China
Room 1001, China Life Tower
16 Chaoyangmenwai Avenue,
Beijing, 100020
Tel.: +86 10 85251111 App. 330
Fax: +86 10 85251001
chinaoffice@tintometer.com
www.lovibond.com
China

Tintometer South East Asia
Unit B-3-12, BBT One Boulevard,
Lebuh Nilam 2, Bandar Bukit Tinggi,
Klang, 41200, Selangor D.E
Tel.: +60 (0)3 3325 2285/6
Fax: +60 (0)3 3325 2287
lovibond.asia@tintometer.com
www.lovibond.com
Malaysia

Tintometer Brazil
Caixa Postal: 271
CEP: 13201-970
Jundiaí – SP
Tel.: +55 (11) 3230-6410
sales@tintometer.com.br
www.lovibond.com.br
Brazil

Tintometer Inc.
6456 Parkland Drive
Sarasota, FL 34243
Tel: 941.756.6410
Fax: 941.727.9654
sales@lovibond.us
www.lovibond.com
USA

Tintometer India Pvt. Ltd.
Door No: 7-2-C-14, 2nd, 3rd & 4th Floor
Sanathnagar Industrial Estate,
Hyderabad, 500018
Telangana
Tel: +91 (0) 40 23883300
Toll Free: 1 800 599 3891/ 3892
indiaoffice@lovibond.in
www.lovibondwater.in
India

Tintometer Spain
Postbox: 24047
08080 Barcelona
Tel.: +34 661 606 770
sales@tintometer.es
www.lovibond.com
Spain

Technical changes without notice
Printed in Germany

Lovibond® and Tintometer® are Trademarks of the Tintometer Group of Companies